

SEA Boards Switch Settings

SW1 - 2D Mono/HVPS				U44 - 2D Grey			
DMA							
1	2	3	4	5	6	7	8
DRQ7	DRQ6	DRQ5	DACK7	DACK6	DACK5	NC	NC

				U45 - 2D Grey			
Interrupt							
1	2	3	4	5	6	7	8
IRQ15	IRQ14	IRQ12	IRQ11	IRQ10	NC	NC	NC

				U10 - 2D Grey			
Minimum Number of Slices							
1	2	3	4	5	6	7	8
1	2	3	4	5	6	7	8

				U17 - 2D Grey			
Blanck Slices at End of Image							
1	2	3	4	5	6	7	8
1	2	3	4	5	6	7	8

SW3 - 2D Mono/HVPS					
Probe Select					
1	2	3	4		
On	Off	On	Off	2D	
Off	On	Off	On	2D2	
Off	On	On	Off	HVPS	
Off	On	Off	On	HVPS	
Off	On	On	Off	2D	MRI

				SW2 - 1D			
Command Select							
1	2	3	4	5	6	7	8
cmd4-	cmd4+	cmd3-	cmd3+	cmd2-	cmd2+	cmd1-	cmd1+
Default: 1-4 off and 5-8 on							

SW1 - 1D			
Aux / Command Select			
1	2	3	4
aux2/cmd4-	aux2/cmd4+	aux1/cmd3-	aux1/cmd3+
Default: all on			